



# Herolab

*Manufacturer of Life Science  
Instruments since 1979*



## *HiGen<sup>®</sup> GT*

**Universal High-Speed Benchtop Centrifuge  
Large Volume up to 4 x 1,000 ml**

### ➤ High-Speed Centrifuges with Flexible Axis!

- This provides the best smooth running.
- The separation of the own centering of the motor and the centering of the installed rotor protects the bearing of the entire drive system.
- Thus, your centrifuge will be maintenance free for a very long time.

Catalog No.	Description	Voltage	Phases	Hz
21 11 150	HiCen GT, Standard for Europe - w./o. refrigeration	230 V	1	50/60 Hz
21 11 151	HiCen GT – w./o. refrigeration	115 V	1	50/60 Hz

- 15 angle rotors, 4 swing-out rotors, max. capacity **4 x 1,000 ml**
- speed: 500 – 14,000 rpm, in steps of 100 rpm
- g-force (RCF): 21,913 x g, in steps of 10 x g
- timer: up to 180 min, in 1 min increments, HOLD for longer runs
- 8 acceleration and deceleration ramps
- maintenance free, brushless induction drive
- stainless steel chamber with additional inner guard ring
- overspeed detection, imbalance detection and cut-off
- motorized lid lock, automatic lid opening, emergency lid lock release
- large color display for both current and preset values of speed/g-force and time



### Additional Options (Available on request for *HiCen® GT* against surcharge)

Cat. No.	Accessories and Options	Dimensions in cm	Dimensions in inches
21 10 675	Safety Castor, Working height with Hicen GT approx. 90 cm (35,43 inches)	(W x D x H) 58 x 69 x 40	(W x D x H) 22,83 x 27,17 x 15,75
20 90 930	Password Protected Start		
20 90 940	GLP Conform Data Management		
20 90 950	Audible Signal in Case of Error		
20 90 970	Operating-Hours Indicator		



### Swing-out Rotors for

*HiCen® GT*

Cat. No.	Type	Max. Capacity	Max. Speed	Max. g-Force
20 97 100	<b>AS 2.6</b>	<b>For 2 x 3 microplates</b> , incl. 2 carriers for plates up to 80 mm high + 2 plate holders	4,000 rpm	1,968 x g
20 97 105	caps	Polysulfone Sealing Caps complete with clips (pack of 2 caps)		
20 90 501	<b>AS 4.500</b>	<b>4 x 500 ml</b> , Basic rotor, buckets to be ordered separately		
20 90 540	carrier	For 4 x 4 microplates, Set of 4 carriers	4,000 rpm	2,719 x g
20 90 502	buckets	4 round buckets for 1 x 500 ml w./o. caps, resp. for adapters	4,500 rpm	3,962 x g
25 35 51	bottles	Bottles 500ml PPCO (98 x 114 mm) with caps, 4 bottles		
20 90 550	Trunnion rings	4 x (7 x 50 ml F), Set of 4 trunnion rings with adapters	4,500 rpm	4,093 x g
20 90 551	Trunnion rings	4 x (14 x 15 ml F), Set of 4 trunnion rings with adapters	4,500 rpm	4,071 x g
20 70 060	<b>AS 4.750</b>	<b>4 x 750 ml</b> , Basic rotor incl. <b>windshield</b> , buckets to be ordered separately		
20 70 400	carrier	For 4 x 4 microplates, Set of 4 carriers	3,700 rpm	2,693 x g
20 70 100	buckets	4 rectangular buckets for tube racks	3,000 rpm	1,962 x g
20 70 300	buckets	4 round buckets for 1 x 750 ml each w./o. caps, resp. for adapters	3,000 rpm	2,133 x g
20 70 050	<b>AS 4.1000</b>	<b>4 x 1000 ml</b> , Basic rotor, buckets to be ordered separately		
20 70 400	carrier	For 4 x 4 microplates, Set of 4 carriers	3,700 rpm	2,693 x g
20 70 100	buckets	4 rectangular buckets for tube racks	2,500 rpm	1,363 x g
20 70 300	buckets	4 round buckets for 1 x 750 ml each w./o. caps, resp. for adapters	2,500 rpm	1,481 x g
20 70 200	buckets	4 round buckets for 1 x 1000 ml each w./o. caps, resp. for adapters	2,500 rpm	1,621 x g
20 70 950	caps	Caps for round buckets 750 ml or 1000 ml, (pack of 2 caps)		
20 70 650	Trunnion rings	4 x (8 x 50 ml F), Set of 4 trunnion rings with adapters	2,500 rpm	1,373 x g
20 70 615	Trunnion rings	4 x (19 x 15 ml F), Set of 4 trunnion rings with adapters	2,500 rpm	1,414 x g

### Angle Rotors (with hermetically sealing lid)

Cat. No.	Type	Max. Capacity	Angle	Max. Speed	Max. g-Force
20 96 100	AF 24.2	24 x 1,5/2,2 ml	45°	14,000 rpm	21,913 x g
20 75 100	AF 72.2	72 x 1,5/2,2 ml	40°	8,000 rpm	9,015 x g
20 94 100	AF 20.10	20 x 10 ml	25°	11,000 rpm	13,257 x g
20 35 00	A 12.24	12 x 14 ml	40°	12,000 rpm	14,972 x g
20 26 100	AF 24.16	24 x 16 ml	28°	9,000 rpm	9,979 x g
20 45 100	AF 12.30	12 x 30 ml	28°	12,000 rpm	17,145 x g
20 93 005	AF 8.50.1 Biosafe	8 x 50 ml	34°	10,000 rpm	12,074 x g
20 93 100	AF 8.50.2	8 x 50 ml	25°	9,000 rpm	8,694 x g
20 98 105	AF 8.50.3 Biosafe	8 x 50 ml Falcon	25°	9,000 rpm	9,418 x g
20 44 100	AF 12.50 Biosafe	12 x 50 ml	34°	7,000 rpm	7,122 x g
20 92 105	AF 6.94 Biosafe	6 x 94 ml	25°	10,000 rpm	10,956 x g
20 71 00	A 8.20	8 x 94 ml	25°	8,000 rpm	8,014 x g
20 22 100	AF 6.250	6 x 250 ml	30°	7,000 rpm	7,943 x g
20 23 100	AF 6.500	6 x 500 ml	30°	7,000 rpm	8,436 x g
20 24 100	AF 4.1000	4 x 1000 ml	22°	4,500 rpm	3,810 x g



For detailed Information about the comprehensive range of [centrifuge tubes und bottles](#) from our production and their possible use in centrifuges of other manufacturers please refer to our catalog "Centrifuge Labware".

**Herolab**

**Tubes and Bottles**

For centrifugation

70 mm wide neck

Ready for your application:

- # from 10 ml up to 1,000 ml Filling capacity
- # from different Plastic materials: PP/CO, PC, PE, PF
- # autoclavable (PP/CO, PC, PF)
- # tested for tightness and max. RC3

© 2014

## Specification

## HiCen® GT

Catalog-No. (230V, 50/60 Hz, 1 Phase)	21 11 150
Maximum Capacity	4 x 1,000 ml
Maximum Speed	14,000 min <sup>-1</sup>
Max. g-Force/RCF (depending on rotor)	21,913 x g
Minimum Speed	> 500 min <sup>-1</sup>
Rotor Detection	APRI™ System = Automatic Positive Rotor Identification for Herolab rotors
Program Memory	30 programs
Display	Color TFT Display for current and preset values
User Interface	Operating panel with buttons and one rotary knob
Preset Time	Up to 180 minutes with increments of 1 min., "HOLD" for longer runs
Preset Speed	From 500 rpm up to max. speed with increments of 100 rpm
Preset g-Force/RCF	With increments of 10 x g
Drive System	Direct self-centering drive system with brushless induction motor
Speed Accuracy	Deviation less than 1 % (typical: 0.5 %)
Acceleration Time (depending on rotor)	30 sec. – 2.5 min.
Deceleration Time (depending on rotor)	40 sec. – 3.5 min.
Acceleration/Deceleration Rates	8 ramps in steps of 30 sec. from 0 to 240 sec. in the range of 0 – 1.000 rpm
Braking System	Dynamic electric brake system
Temperature Set Range	n.a.
Cooling at Standstill	n.a.
Temperature Control Accuracy	n.a.
Temperature Control System	n.a.
Sample Protection	n.a.
Overtemperature Protection	n.a.
Refrigeration System	n.a.



Safety Features	<ul style="list-style-type: none"> <li>Overspeed detection</li> <li>Imbalance detection and cut-off</li> <li>Four-point lid fixing</li> <li>Automatic motor driven lid lock with emergency lid lock release</li> <li>Stainless steel chamber with additional inner guard ring</li> </ul>
-----------------	--

Dimensions (W x D x H): in cm (in inches)	57.5 x 76.0 x 50.0 (22.6 x 29.9 x 19.7)
--	--

Shipping Dimensions in cm (in inches)	105 x 95 x 60 (41,3 x 37,4 x 23.6)
--	---------------------------------------

Net Weight	112 kg (247 lbs)
------------	------------------

Gross Weight	132 kg (291 lbs)
--------------	------------------



WEEE-Reg. No.  
DE66734561



All products built by Herolab carry the CE sign.  
Herolab has its own quality management system according to ISO 9001.

# Herolab




Herolab GmbH Laborgeräte

Ludwig-Wagner-Str. 12  
69168 Wiesloch  
GERMANY  
Tel.: +49 (0) 6222 – 5802-0  
Fax: +49 (0) 6222 – 5802-34  
E-Mail: info@herolab.de  
Internet: www.herolab.de



## Tubes and Adapters for Rotors of *HiGen® GT*

for Rotors **AF 24.2 + AF 72.2**

Tubes	2.0-2.2 ml	1.5 ml	0.5-0.7 ml	0.25-0.4 ml	0.2 ml
Tubes Ø x Length (mm)	11/13 x 40	11/13 x 40	8/10 x 28-31	Ø 6 mm	PCR® Tubes
Material of Tubes + Herolab Cat. No.	Tubes not available from Herolab	Tubes not available from Herolab	Tubes not available from Herolab	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter	---	---			
Herolab Cat. No.	---	---	25 50 06	25 50 05	25 50 07
Tubes per Adapter / Rotor	- / 24 resp. 72	- / 24 resp. 72	1 / 24 resp. 72	1 / 24 resp. 72	1 / 24 resp. 72



Cat. No. 20 75 100



Cat. No. 20 96 100



Cat. No. 25 24 01 / 11 / 41 / 51

Cat. No. 20 94 100



Cat. No. 25 25 70 / 25 80 / 26 60 / 25 90 / 26 90

for Rotor




### AF 20.10

### A 12.24

Tubes	10 ml AOR	13.5 ml
Tubes Ø x Length (mm)	16 x 80	16 x 76
Material of Tubes + Herolab Cat. No.	PC: 25 24 01 PPCO: 25 24 11 PE: 25 24 41 PF: 25 24 51	PPCO: 25 24 70
Necessary Adapter	---	---
Herolab Cat. No.	---	---
Tubes per Adapter / Rotor	- / 20	- / 20



Cat. No. 20 35 00

14 ml	5 ml Glass	4.5 / 5 ml	1.5 ml
18 x 75	12 x 75	12 x 75	11/13 x 40
PC: 25 25 70* PE: 25 26 60* PF: 25 25 90* PPCO: 25 26 90*	Tubes not available from Herolab	PP: 25 22 30	Tubes not available from Herolab
---			
---	25 56 40	25 50 73	25 50 71
- / 12	1 / 12	1 / 12	1 / 12

Cat. No. 25 27 00 / 20 / 40 / 51



Cat. No. 20 26 100



\* Snap-on caps for 14 ml and 16 ml tubes available (Cat No. 25 45 50 / 25 45 51)

Cat. No. 25 29 51 / 61 / 65 / 71



for Rotor

### AF 24.16

### AF 12.30

Tubes	16 ml with rim	12/15 ml Glass	10 ml Glass	4.5 / 5 ml	3 ml Glass	1.5 ml
Tubes Ø x Length (mm)	18 x 100	18 x 102	13 x 100	12 x 75	10 x 75	11/13 x 40
Material of Tubes + Herolab Cat. No.	PC: 25 27 00* PPCO: 25 27 20* PE: 25 27 40* PF: 25 27 51*	15 ml HS: 25 10 91	Tubes not available from Herolab	PP: 25 22 30	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter	---					
Herolab Cat. No.	---	25 58 20	25 56 00	25 50 72	25 55 80	25 50 70
Tubes per Adapter / Rotor	- / 24	1 / 24	1 / 24	1 / 24	1 / 24	1 / 24










Cat.-or. 25 30 20 + 25 46 85

30 ml AOR
25 x 92
PC: 25 29 51 PPCO: 25 29 61 PE: 25 29 65 PF: 25 29 71
---
---
- / 12

## Tubes and Adapters for *HiGen<sup>®</sup> GT*

for Rotor **AF 8.50.1 + AF 8.50.2 + A 12.17**

Tubes	50ml AOR	50 ml with rim	50 ml with cap	30 ml AOR	16 ml with rim	12/15 ml Glass	10ml AOR	4.5/5 ml
Tubes Ø x Length (mm)	29 x 103	29 x 103	29 x 103	25 x 92	18 x 100	18 x 102	16 x 80	12 x 75
Material of Tubes + Herolab Cat. No.	PC: 25 32 01 PPCO: 25 32 11 PE: 25 32 15 PF: 25 32 21	PC: 25 31 00 PPCO: 25 31 60 PE: 25 31 30 PF: 25 31 51	PC: 25 31 05 PPCO: 25 31 45	PC: 25 29 51 PPCO: 25 29 61 PE: 25 29 65 PF: 25 29 71	PC: 25 27 00 PPCO: 25 27 20 PE: 25 27 40 PF: 25 27 51	Tubes not available from Herolab	PC: 25 24 01 PPCO: 25 24 11 PE: 25 24 41 PF: 25 24 51	PP: 25 22 30
Necessary Adapter	---	---	---					
Herolab Cat. No.	---	---	---	25 50 61	25 50 50	25 55 50	25 50 30	25 50 11
Tubes per Adapter / Rotor	- / 8 resp. 12	- / 8 resp. 12	- / 8 resp. 12	1 / 8 resp. 12	1 / 8 resp. 12	1 / 8 resp. 12	1 / 8 resp. 12	2 / 16 resp. 24

Tubes	1.5 ml	1 ml
Tubes Ø x Length (mm)	11 x 40	7 x 50
Material of Tubes + Herolab Cat. No.	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter		
Herolab Cat. No.	25 50 10	25 50 00
Tubes per Adapter / Rotor	2 / 16 resp. 24	4 / 32 resp. 48





Cat. No. 25 32 01 / 11 / 15 / 21

Cat. No. 25 31 05 / 45

Cat. No. 25 31 00 / 60 / 30 / 51

for Rotor **AF 8.50.3**

Tubes	50 ml Falcon	30 ml AOR	15 ml Falcon
Tubes Ø x Length (mm)	29 x 118	25 x 92	17 x 122
Material of Tubes + Herolab Cat. No.	Tubes not available from Herolab	PC: 25 29 51 PPCO: 25 29 61 PE: 25 29 65 PF: 25 29 71	Tubes not available from Herolab
Necessary Adapter	---		
Herolab Cat. No.	---	25 50 62	25 53 80
Tubes per Adapter / Rotor	- / 8	1 / 8	1 / 8



Cat. No. 20 44 00



Cat. No. 20 93 000



Kat.-Nr. 20 9§ 100




Cat. No. 20 98 100

## Tubes and Adapters for *HiGen® GT*

### for Rotor **AF 6.250**

Tubes/Bottles	250 ml	250 ml	200ml	100/120 ml	30 ml Glass	16 ml with rim	12/15 ml	10 ml
Tubes Ø x Length (mm)	62 x 122	62 x 122	60 x 115	40 x 115	24 x 105	18 x 100	18 x 102	13 x 100
Material of Tubes + Herolab Cat. No.	PC: 25 34 01 PPCO: 25 34 21 PE: 25 34 31 PF: 25 34 39 PP black: 25 34 12	round bottom: PC: 25 34 41 PPCO: 25 34 43 PPCO: 25 34 45 PE: 25 34 47 PF: 25 34 49	Tubes not available from Herolab	100 ml: PC: 25 33 00	Tubes not available from Herolab	PC: 25 27 00 PPCO: 25 27 20 PE: 25 27 40 PF: 25 27 51	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter	---							
Herolab Cat. No.	---	25 51 90	25 58 10	25 51 70	25 51 60	25 51 30	25 51 30	25 51 10
Tubes per Adapter / Rotor	-	1 / 6	1 / 6	1 / 6	3 / 18	5 / 30	5 / 30	9 / 54

Tubes	4.5/5 ml	3 ml	1.5 ml
Tubes Ø x Length (mm)	12 x 75	10 x 75	11/13 x 40
Material of Tubes + Herolab Cat. No.	PP: 25 22 30	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter			
Herolab Cat. No.	25 51 00	25 50 90	25 52 10
Tubes per Adapter / Rotor	9 / 54	15 / 90	8 / 48



Cat. No. 25 34 01 / 21 / 31 / 39 / 12

Cat. No. 25 34 41 / 43 / 45 / 47 / 49



Cat. No. 20 22 100

### for Rotor **AF 6.500**

Bottles	500 ml
Ø x Length (mm)	70 x 165
Material of Tubes + Herolab Cat. No.	PC: 25 35 01* PPCO: 25 35 21* PE: 25 35 31* PP black: 25 35 12*
Necessary Adapter	---
Herolab Cat. No.	- / 6
Tubes per Adapter / Rotor	



\* For safety reasons always use 500 ml bottles with collar support ring. (Catalog No. 25 52 41)!



Cat. No. 25 35 01 / 21 / 31 / 12

### **AF 4.1000**

1000 ml
(100 x 181)
(PPCO: 25 37 50** for rotor 20 24 110 only)
98 x 175
Cap with Support Ring: PC: 25 35 65** PPCO: 25 35 75**
98 x 190
PPCO: 25 37 80**
---
- / 4



Cat. No. 25 35 65 / 75

Cat. No. 25 37 80

Cat. No. 25 37 50

#### Widemouth Bottle

\*\* Use the Ø 98 mm bottles in fixed angle rotors for 98 mm Ø bottles only with cover including support ring. For rotors for 100 mm Ø bottles the stainless-steel cup 25 13 511 should be used as an adapter. Attention: Then pay attention to the total weight per bore! The 100 mm bottles are wide-mouth bottles with an opening of 70 mm.



Cat. No. 20 24 100



## Tubes and Adapters for *HiGen® GT*

for Rotor **AF 6.94 + A 8.20**

Cat. No. 20 71 00



Cat. No. 20 92 100

Cat. No. 25 32 42 / 45 / 47 / 49



Cat. No. 25 32 90 + 25 46 82

94 ml	80 ml AOR
<b>38 x 102</b>	38 x 105
thick-walled PC: 25 32 80 PPCO: 25 32 60 thick-w. w. thread PC: 25 32 90	PC: 25 32 41 PPCO: 25 32 45 PE: 25 32 47 PF: 25 32 49
---	---
---	---
<b>- / 6 bzw. 8</b>	<b>- / 6 bzw. 8</b>

Adapter for tubes (38.5 ml), 17 ml, 13.5 ml, 6.5 ml and 4.4 ml are available on request.



Cat. No. 25 30 20 + 25 46 85

for Rotor **AS 2.6**

Plates	Microtiter Plates	Deepwell Plates
Total Height of Plates	up to 80 mm	up to 80 mm
Material of Tubes + Herolab Cat. No.	Plates not available from Herolab	Plates not available from Herolab
Necessary Adapter	---	---
Herolab Cat. No.	---	---
Tubes per Adapter / Rotor	- / 2-6	- / 2-4


**AS 4.750 + AS 4.1000**  
with set of Microtiter Plate carriers

Microtiter Plates	Deepwell Plates
up to 64 mm	up to 64 mm
Plates not available from Herolab	Plates not available from Herolab
---	---
---	---
- / 4-12	- / 4-8
---	---



Cat. No. 20 70 400

Cat. No. 20 97 100

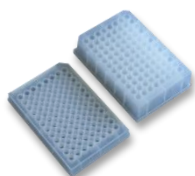

Cat. No. 20 97 105



Cat. No. 20 70 060





Cat. No. 20 70 050





## Tubes and Adapters for *HiCen<sup>®</sup>GT*

for Rotor **AS 4.500**

Tubes/Bottles	500 ml	50 ml Falcon	15 ml Falcon
Tubes Ø x Length (mm)	98 x 114	29 x 118	17 x 122
Material of Tubes + Herolab Cat. No.	PPCO: 25 35 51	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter	---		
Herolab Cat. No.	---	<b>20 90 530</b>	<b>20 90 535</b>
Tubes per Adapter / Rotor	20 70 060 / 4	5 / 20	8 / 32
(Pack of 2 Adapters)	---		

Cat. No. 20 90 501 + 20 90 511  
+ 20 90 530 + 20 90 535



Cat. No. 25 35 51

for Rotor **AS 4.750 + AS 4.1000**  
with set of rectangular buckets for tube racks

Tubes/Bottles	250 ml	100 ml	50 ml Falcon	50ml AOR	50 ml with rim	50 ml with cap	30 ml AOR	15 ml Falcon
Tubes Ø x Length (mm)	62 x 122	40 x 115	29 x 118	29 x 103	29 x 103	29 x 103	25 x 92	17 x 122
Material of Tubes + Herolab Cat. No.	PC: 25 34 01 PPCO: 25 34 21 PE: 25 34 31 PF: 25 34 39 PPblack: 25 34 12	PC: 25 33 00	Tubes not available from Herolab	PC: 25 32 01 PPCO: 25 32 11 PE: 25 32 15 PF: 25 32 21	PC: 25 31 00 PPCO: 25 31 60 PE: 25 31 30 PF: 25 31 51	PC: 25 31 05 PPCO: 25 31 45	PC: 25 29 51 PPCO: 25 29 61 PE: 25 29 65 PF: 25 29 71	Tubes not available from Herolab
Necessary Adapter								
Herolab Cat. No.	<b>20 70 110</b>	<b>20 70 115</b>	<b>20 70 130</b>	<b>20 70 125</b>			<b>20 70 135</b>	<b>20 70 145</b>
Tubes per Adapter / Rotor	4	4	2 / 8	2 / 8			3 / 12	6 / 24

Cat. No. 20 70 100

Kat.-Nr. 25 34 01 / 21 / 31 / 39 / 12



Cat. No.  
25 32 01 / 11  
25 32 15 / 21

Cat. No.  
25 31 00 / 60  
25 31 30 / 51



Cat. No.  
25 31 05 / 45

Cat. No.  
25 29 51 / 61  
25 32 65 / 71

## Tubes and Adapters for *HiGen<sup>®</sup> GT*



Kat.-Nr. 25 35 62 / 72      Kat.-Nr. 25 35 80 / 81



Kat.-Nr. 20 70 200      Kat.-Nr. 20 70 300

for Rotor **AS 4.750 + AS 4.1000**  
with set of round buckets for bottles or adapters

Tubes/Bottles	1000 ml	750 ml	500 ml	250 ml	250 ml	100 ml	50 ml Falcon
Tubes Ø x Length (mm)	98 x 175	98 x 152	70 x 165	62 x 122	Ø 63	40 x 115	29 x 118
Material of Tubes + Herolab Cat. No.	PC: 25 35 62 PPCO: 25 35 72	PC: 25 35 80 PPCO: 25 35 81	PC: 25 35 01* PPCO: 25 35 21* PE: 25 35 31* PP black: 25 35 12*	PC: 25 34 01 PPCO: 25 34 21 PE: 25 34 31 PF: 25 34 39 PP black: 25 34 12	Conical bottles  Bottles not available from Herolab	PC: 25 33 00	Tubes not available from Herolab
Necessary Adapter	--- Can only be used in 1000 ml buckets No. 20 70 200 of rotor AS 4.1000 directly	--- Can only be used in 750 ml buckets No. 20 70 300 of rotor AS 4.1000 directly					
Herolab Cat. No.	---	---	20 70 210	20 70 220	20 70 215	20 70 225	20 70 230
Tubes per Adapter / Rotor	- / 4	- / 4	1 / 4	1 / 4	1 / 4	3 / 12	5 / 20

\* For safety reasons always use 500 ml bottles with cover support ring (Cat. No. 25 52 41)!



Tubes	50ml AOR	50 ml with rim	50 ml with cap	30 ml AOR	15ml	15 ml Falcon	10ml AOR
Tubes Ø x Length (mm)	29 x 103	29 x 103	29 x 103	25 x 92	17 x 87/119	17 x 122	16 x 80
Material of Tubes + Herolab Cat. No.	PC: 25 32 01 PPCO: 25 32 11 PE: 25 32 15 PF: 25 32 21	PC: 25 31 00 PPCO: 25 31 60 PE: 25 31 30 PF: 25 31 51	PC: 25 31 05 PPCO: 25 31 45	PC: 25 29 51 PPCO: 25 29 61 PE: 25 29 65 PF: 25 29 71	Tubes not available from Herolab	Tubes not available from Herolab	PC: 25 24 01 PPCO: 25 24 11 PE: 25 24 41 PF: 25 24 51
Necessary Adapter							
Herolab Cat. No.	20 70 235			20 70 240	20 70 250	20 70 245	20 70 255
Tubes per Adapter / Rotor	7 / 28			8 / 32	18 / 72	12 / 48	20 / 80

Tubes	7 ml	7 ml	5 ml	0.5 ml
Tubes Ø x Length (mm)	Ø 13.5	Ø 16.5	Ø 13.5	Ø 9.5
Material of Tubes + Herolab Cat. No.	Tubes not available from Herolab	Tubes not available from Herolab	Tubes not available from Herolab	Tubes not available from Herolab
Necessary Adapter				
Herolab Cat. No.	20 70 260	20 70 265	20 70 270	20 70 275
Tubes per Adapter / Rotor	32 / 128	21 / 84	35 / 140	60 / 240

Cat. No. 25 35 01 / 21 / 31 / 12      Cat. No. 25 34 01 / 21 / 31 / 39 / 12



Cat. No. 25 32 01 / 11 / 15 / 21 + 25 31 05 / 45 / 00 / 60 / 30 / 51



Cat. No. 25 32 01 / 11  
25 32 15 / 21



Cat. No. 25 29 51 / 61  
25 32 65 / 71

Herolab Tubes for Rotors of *HiGen<sup>®</sup> GT*

Cat. No.	Description	Ø x Length (mm)	Pieces/Pack
25 22 30	Tubes 4.5 ml PP-RB	12 x 75	1000
25 24 01	Tubes 10 ml PC-AOR with screw cap	16 x 80	10
25 24 11	Tubes 10 ml PPCO-AOR with screw cap	16 x 80	10
25 24 41	Tubes 10 ml PE-AOR with screw cap	16 x 80	10
25 24 51	Tubes 10 ml PF-AOR with screw cap	16 x 80	2
25 24 70	Tubes 13.5 ml PPCO-RB	16 x 76	10
25 25 70	Tubes 14 ml PC-RB	18 x 75	10
25 25 80	Tubes 14 ml PPCO-RB	18 x 75	10
25 25 90	Tubes 14 ml PF-RB	18 x 75	2
25 26 60	Tubes 14 ml PE-RB	18 x 75	10
25 26 90	Tubes 14 ml PPCO-RB	18 x 75	10
25 27 00	Tubes 16 ml PC-RB/FL	18 x 100	10
25 27 20	Tubes 16 ml PPCO-RB/FL	18 x 100	10
25 27 40	Tubes 16 ml PE-RB/FL	18 x 100	10
25 27 51	Tubes 16 ml PF-RB/FL	18 x 100	2
25 29 51	Tubes 30 ml PC-AOR with screw cap	25 x 92	10
25 29 61	Tubes 30 ml PPCO-AOR with screw cap	25 x 92	10
25 29 65	Tubes 30 ml PE-AOR with screw cap	25 x 92	10
25 29 71	Tubes 30 ml PF-AOR with screw cap	25 x 92	2
25 30 20	Tubes 38.5 ml PC thick-walled with thread UC	25 x 89	50
25 30 40	Tubes 38.5 ml PC thick-walled UC	25 x 89	50
25 30 50	Tubes 38.5 ml PPCO thin-walled UC	25 x 89	50
25 30 60	Tubes 38.5 ml PPCO thick-walled UC	25 x 89	50
25 31 00	Tubes 50 ml PC-RB/FL	29 x 103	10
25 31 05	Tubes 50 ml PC-RB/FL with Loop Cap	29 x 103	10
25 31 30	Tubes 50 ml PE-RB/FL	29 x 103	10
25 31 45	Tubes 50 ml PPCO-RB/FL with Loop Cap	29 x 103	10
25 31 51	Tubes 50 ml PF-RB/FL	29 x 103	2
25 31 60	Tubes 50 ml PPCO-RB/FL	29 x 103	10
25 32 01	Tubes 50 ml PC-AOR with screw cap	29 x 103	10
25 32 11	Tubes 50 ml PPCO-AOR with screw cap	29 x 103	10
25 32 15	Tubes 50 ml PE-AOR with screw cap	29 x 103	10
25 32 21	Tubes 50 ml PF-AOR with screw cap	29 x 103	2
25 32 41	Tubes 80 ml PC-AOR with screw cap	38 x 105	10
25 32 45	Tubes 80 ml PPCO-AOR with screw cap	38 x 105	10
25 32 47	Tubes 80 ml PEC-AOR with screw cap	38 x 105	10
25 32 49	Tubes 80 ml PF-AOR with screw cap	38 x 105	2
25 32 50	Tubes 94 ml PPCO-RB	38 x 106	10
25 32 60	Tubes 94 ml PPCO thick-walled UC	38 x 102	25
25 32 80	Tubes 94 ml PC thick-walled UC	38 x 102	25
25 32 90	Tubes 94 ml PC thick-walled with thread UC	38 x 102	25
25 33 00	Tubes 100 ml PC-RB/G	40 x 115	2
25 33 10	Tubes 120 ml PP-RB	40 x 115	10
25 34 01	Bottles 250 ml PC-FB with screw cap	62 x 122	6
25 34 12	Bottles 250 ml PPCO black FB with screw cap	62 x 122	6
25 34 21	Bottles 250 ml PPCO-FB with screw cap	62 x 122	6
25 34 31	Bottles 250 ml PE-FB with screw cap	62 x 122	6
25 34 39	Bottles 250 ml PF-FB with screw cap	62 x 122	2
25 34 41	Bottles 250 ml PC-RB with screw cap	62 x 122	6
25 34 43	Bottles 250 ml PPCO-RB with screw cap	62 x 122	6
25 34 45	Bottles 250 ml PPCO-RB with screw cap	62 x 122	6
25 34 47	Bottles 250 ml PE-RB with screw cap	62 x 122	6
25 34 49	Bottles 250 ml PF-RB with screw cap	62 x 122	2
25 34 81	Bottles 290 ml PC-FB with screw cap	62 x 137	6
25 34 87	Bottles 290 ml PE-FB with screw cap	62 x 137	6
25 35 01	Bottles 500 ml PC-FB with screw cap	70 x 165	6

Cat. No.	Description	Ø x Length (mm)	Pieces/Pack
25 35 12	Bottles 500 ml PPCO black-FB with screw cap	70 x 165	2
25 35 21	Bottles 500 ml PPCO-FB with screw cap	70 x 165	6
25 35 31	Bottles 500 ml PE-FB with screw cap	70 x 165	6
25 35 51	Bottles 500 ml PPCO-FB with screw cap (for AS 4.500 only)	98 x 114	4
25 35 65	Bottles 1000 ml PC-FB with screw cap with neck support ring	98 x 175	4
25 35 75	Bottles 1000 ml PPCO-FB with screw cap with neck support ring	98 x 175	4
25 37 50	Bottles 1000 ml PPCO-FB with sealing cap	100 x 181	4
25 37 80	Bottles 1000 ml PPCO-FB with sealing cap	98 x 190	4
25 45 50	Snap-on Caps for Tubes 14 ml and 16 ml (except for PE)		10
25 45 51	Snap-on Caps for Tubes 14 ml and 16 ml PE		10
25 46 82	Screw cap, aluminum, for UC tubes 94 ml		1
25 46 85	Screw cap, aluminum, for UC tubes 38,5 ml		1
25 52 41	Support ring for bottles 500 ml		6



UC	Ultra Centrifuge Tubes	PPCO	Polypropylene Copolymer
AOR	Advanced Oak Ridge	CB	Conical Bottom
PC	Polycarbonate	FB	Flat Bottom
PE	Polyethylene	RB	Round Bottom
PF	Polyflor®	FL	Flange (Rim)

Herolab reserves the right to change technical specification and prices without prior notice.